


**ACIDE PERACETIQUE 0-50PPM BTE 100 BANDELETTES**

Référence

0299004

Low Level Peracetic Acid Test Strip, 50ppm

1. Remove one test strip from the vial, being careful not to touch the test pad with your fingers.
2. Gently swirl test strip in small sample of solution being tested for 2 seconds.
3. Remove strip with pad face up. Do not shake off excess sample.
4. Compare the test pad to the color chart after 10 seconds.

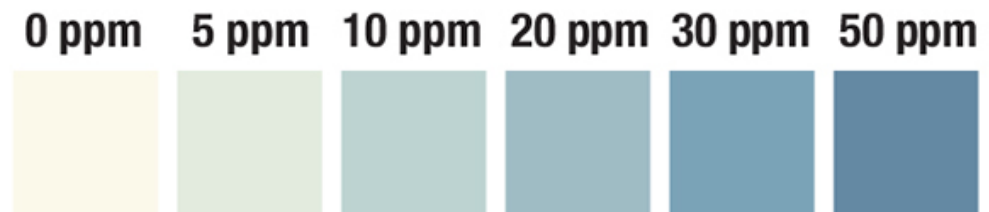

*Ces spécifications ne peuvent être considérées que comme des informations générales et ne dispenseront pas l'utilisateur de son obligation de s'assurer que le produit est adapté pour l'utilisation prévue. En cas de doute, veuillez nous contacter.*

FT 0299004

avril 24