BAIRD PARKER AGAR + NEUTRALIZING

Selective medium for *Staphylococcus aureus* isolation with inactivation of disinfectants.

TYPICAL FORMULA (g/l)

Tryptone	10.0
Beef Extract	5.0
Yeast Extract	1.0
Glycine	12.0
Sodium Pyruvate	10.0
Lithium Chloride	5.0
Agar	17.0
Sodium Phosphate Bibasic	0.4
Lecithin	0.03
L-Histidine	0.01
Sodium Thiosulfate	0.08
Tween 80	0.3
Citric Acid	0.1
Egg Yolk	35.0 ml
Potassium Tellurite	0.1
Final pH 7.0 ± 0.2	

DESCRIPTION

BAIRD PARKER AGAR + NEUTRALIZING is a selective medium for *Staphylococcus aureus* isolation with inactivation of disinfectants.

PRINCIPLE

Tryptone and beef extract are a most satisfactory source of nitrogen and contain a mix of peptides and free amino acids. Yeast extract is a source of amino acids and vitamin B. Sodium pyruvate is incorporated in order to stimulate the growth of *S. aureus* without destroying the selectivity. The tellurite is toxic to egg yolk-clearing strains other than *S. aureus* and imparts a black color to the colonies of *S. aureus*. The egg yolk, in addition to be an enrichment, aids in the identification process by demonstrating lecithinase activity. Glycine and lithium chloride have inhibitory action for organisms other than *S. aureus*. Lecithin and Tween 80 are the ingredients of the neutralizing which eliminates the bactericidal activity of ammonium quaternary compounds contained in sanitizers.

TECHNIQUE

- Remove the first wrapper of the packet and the inner one.
- Remove the cap of the plate and gently push the agar surface on a level surface to examine for at least 10 seconds. Otherwise streak the sampling swab onto the surface of the medium.
- Cover the plate with the cap.
- Incubate for 24-48 hours at 36±1°C.

INTERPRETATION of RESULTS

Observe for the formation of Units Forming Colonies (UFC).

STORAGE

2-8°C away from light, until the expiry date on the label or until signs of deterioration or contamination are evident.

WARNING and PRECAUTIONS

The product is not classified as hazardous by current legislation and does not contain harmful substances in concentrations of ≥1%. The product is designed for *In vitro* diagnostic use and must be used only by properly trained operators.

DISPOSAL of WASTE

Disposal of waste must be carried out according to national and local regulations in force.

REFERENCES

1. Baird Parker, A.C.. (1962). *An. J. Appl. Bacteriol.* 25:12-19.
2. Baird Parker, A.C. (1969) *Isolation methods for microbiologists.* Shapton,. & Gould ed. London: Academic Press.
3. ISO 6888-1:1999. *Part 1: Technique using Baird-Parker agar medium.*

Liofilchem s.r.l.

Via Scozia, Zona industriale - 64026 Roseto degli Abruzzi (Te) Italy - Tel. +39.0858930745 - Fax +39.0858930330
Web site: <http://www.liofilchem.net> - E-mail: liofilchem@liofilchem.net

PRODUCT SPECIFICATIONS

NAME

BAIRD PARKER AGAR + NEUTRALIZING

PRESENTATION

Rodac plates containing 17+/-1 ml of medium.

STORAGE

2-8°C

PACKAGING

Code	Content	Packaging
15331	20 plates	1 plate in blister 20 plates in cardboard box 1 instructions sheet - Outside label

pH OF THE MEDIUM

7.0 ± 0.2

USE

BAIRD PARKER AGAR + NEUTRALIZING is a selective medium for *Staphylococcus aureus* isolation with inactivation of disinfectants.

TECHNIQUE

Refer to technical sheet of the product.

APPEARANCE of the MEDIUM

Yellow medium, opalescent.

SHELF LIFE

4 months

QUALITY CONTROL

- Control of general characteristics, label and print
- Sterility control
7 days at 25 ± 1°C, in aerobiosis
7 days at 36 ± 1°C, in aerobiosis
- Microbiological control
Inoculum for productivity: 10-100 UFC/ml
Inoculum for selectivity: 10⁴-10⁵ UFC/ml
Inoculum for specificity: ≤ 10⁴ UFC/ml
Incubation conditions: 48 h at 36 ± 1°C, in aerobiosis

Microorganism		Growth	Characteristics	Lecithinase production (halos)
<i>Escherichia coli</i>	ATCC 25922	Inhibited		-
<i>Bacillus subtilis</i>	ATCC 6633	Poor	Brown colonies	-
<i>Proteus mirabilis</i>	ATCC 25933	Good	Brown colonies	-
<i>Staphylococcus aureus</i>	ATCC 25923	Good	Black colonies	+
<i>Staphylococcus epidermidis</i>	ATCC 14990	Poor	Black colonies	-

TABLE OF SYMBOLS

LOT Batch code	 Caution, consult accompanying documents	 Manufacturer	 Contains sufficient for <n> tests
REF Catalogue number	 Temperature limitation	 Use by	

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