## Acetamide Broth

Liquid medium for detection of *P. aeruginosa* in water samples, according to ISO 16266.

### DESCRIPTION

Acetamide Broth is a liquid medium used for confirmation of *Pseudomonas aeruginosa* with the membrane filtration technique.

This medium complies with ISO 16266 for examination of water with low background flora such as waters intended for human consumption or pool waters.

| TYPICAL FORMULA                | (g/l)  |
|--------------------------------|--------|
| Potassium Di-hydrogenphosphate | 1.0    |
| Magnesium Sulfate, anhydrous   | 0.2    |
| Acetamide                      | 2.0    |
| Sodium Chloride                | 0.2    |
| Sodium Molybdate               | 0.005  |
| Iron Sulfate, heptahydrate     | 0.0005 |
| Final pH 7.0 ± 0.5 at 25°C     |        |

### METHOD PRINCIPLE

Potassium phosphate is the buffering agent. Magnesium sulfate, sodium molybdate and iron sulfate are inorganic salts essential for organisms growth. Acetamide is the sole source of carbon in the medium. Sodium chloride maintains the osmotic balance of the medium. Acetamide is deaminated by gram-negative non-fermenting bacteria to produce ammonia.

### PREPARATION

Dehydrated medium Suspend 3.4 g of the powder in 1 liter of distilled or deionized water. Mix well. Heat to boil shaking frequently until completely dissolved. Dispense into tubes or other final containers. Sterilize in autoclave at 121°C for 15 minutes.

### TEST PROCEDURE

Filter volumes of diluted or undiluted water sample through 0.45 µm sterile membranes. Place each membrane on a plate of *Pseudomonas* CN Agar (ref. 163592) and incubate at 36 ± 2°C for 40-48 hours. Count all colonies that produce blue/green (pyocyanin) color as confirmed *P. aeruginosa*.

Acetamide Broth is used for confirming fluorescent non-pyocyanin producing colonies and reddish brown pigmented colonies:

- Transfer all colonies to Nutrient Agar (ref. 10044) and incubate at 36 ± 2°C for 20-24 h;
- Inoculate a tube of Acetamide Broth with the subculture and incubate at 36 ± 2°C for 20-24 h;
- Add 1-2 drops of Nessler's Reagent (ref. 80284) and examine for production of ammonia.

### INTERPRETING RESULTS

The formation of a yellow to brick red color indicates ammonia production. A negative reaction shows no color change.

Other confirmatory tests, such as Oxidase Test (ref. 88029, 88003 or 88004) and King's B Medium (ref. 11072), may be required.

### APPEARANCE

Dehydrated medium: free-flowing, homogeneous, white to light beige.

Prepared medium: Colourless, limpid solution.

### STORAGE

The powder is very hygroscopic, store the powder at 10-30°C, in a dry environment, in its original container tightly closed. Store tubes at 2-8°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

### SHELF LIFE

Dehydrated medium: 4 years.

Medium in tubes: 1 year.

**QUALITY CONTROL**

The medium is inoculated with the microbial strains indicated in the QC table.

Inoculum for productivity:  $\leq 100$  CFU.

Incubation conditions: aerobically at  $36 \pm 2^\circ\text{C}$  for 20-24 hours.

**QC Table.**

| Microorganism                 |             | Growth | Acetamide Test |
|-------------------------------|-------------|--------|----------------|
| <i>Pseudomonas aeruginosa</i> | ATCC® 27853 | Good   | Positive       |
| <i>Pseudomonas aeruginosa</i> | ATCC® 10145 | Good   | Positive       |
| <i>Pseudomonas aeruginosa</i> | ATCC® 9027  | Good   | Positive       |
| <i>Escherichia coli</i>       | ATCC® 25922 | Good   | Negative       |

**WARNING AND PRECAUTIONS**

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use and must be used only by properly trained operators.

**DISPOSAL OF WASTE**

Disposal of waste must be carried out according to national and local regulations in force.

**BIBLIOGRAPHY**

1. EN ISO 16266: 2008. Water quality – Detection and enumeration of *Pseudomonas aeruginosa* – Method by membrane filtration (ISO 16266:2006).

| PRESENTATION    | Format            | Packaging       | Ref.   |
|-----------------|-------------------|-----------------|--------|
| Acetamide Broth | Tube (20x100mm)   | 20 x 5 ml tubes | 24154  |
| Acetamide Broth | Dehydrated medium | 500 g of powder | 610392 |

**TABLE OF SYMBOLS**

|                             |                                                                                                             |                                                                                                                       |                                                                                                                          |                                                                                                                 |
|-----------------------------|-------------------------------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------|
| <b>LOT</b> Batch code       |  Keep away from sunlight |  Manufacturer                      |  Use by                               |  Fragile, handle with care |
| <b>REF</b> Catalogue number |  Temperature limitation  |  Contains sufficient for <n> tests |  Caution, consult Instruction For Use |  Do not reuse              |

This document is available from the online Support Center:

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