



IPA Surface Disinfectant Wipes



3 products in this range

Powerful, fast-acting disinfection.

- Broad-spectrum efficacy - helping you maintain your environment with confidence
- Rapid action without leaving any visible residue behind
- 70% isopropyl alcohol (IPA), the perfect choice for water-sensitive areas
- EU BPR Union approved



IPA Surface Disinfectant Wipes



Key Highlights

- For use disinfecting food areas
- Quat & PHMB Free
- HALAL certified
- Suitable for use within dry food processing and manufacturing
- Blue, visually detectable wipe material
- EU Union Authorisation (PT2, PT4) - EU-0020463-0000
- 2l canister - 27% plastic reduction compared with previous version
- Tritex material, canister, buckets, lids, and labels are fully recyclable
- Cardboard used in outer packaging is FSC approved

How to use

1



Risk assessment

Please follow your agreed risk assessment policy guidelines regarding the use of PPE.


4



S-shaped motion

Wipe in an S-shaped motion from clean to dirty, applying even pressure with the wipe flat. Do not rewipe the same area. Discard the wipe if it becomes soiled or dry, and take a new one from the pack.

2

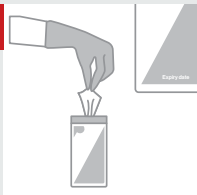

Wipe selection

Ensure the wipes are still within the expiry date before use. Close the pack after selecting your wipes to prevent contamination.

5



Discard waste

Discard used wipes in the designated waste bin, following your local guidelines.


3



Compatibility

Ensure you use the correct wipe for the surface or equipment, following protocols and compatibility guidelines.

6

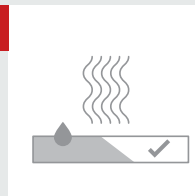

Air dry

Make sure the surface stays wet for the required contact time as per the packaging instructions. Ensure the surface is fully dry before use.

Technical details

Chemical Formulation

Paltx IPA Surface Disinfectant Wipes contain 70% Isopropyl Alcohol (IPA).

A 70% alcohol solution is optimal for disinfection, the added water slows evaporation, increases surface contact time, and helps the alcohol penetrate microbial cell membranes.

Substrate*

Pal's proprietary Tritex® material is made by bonding layers of spunbond and meltblown polypropylene fibres together in a high heat process.

The addition of a hydrophilic coating means optimum and consistent liquid discharge.

Tritex® is 100% synthetic.

*Excludes W84230T.

Substrate key features



High Strength



Excellent solution retention



Very low linting



Efficient solution release

Efficacy details

Standard	Organism Group	Organism(s)	Contact Time	Conditions
EN 1276	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i> <i>Escherichia coli</i>	1 minute	Clean
EN 1276	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i> <i>Escherichia coli</i>	1 minute	Dirty
EN 13727	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i>	1 minute	Dirty
EN 13727	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i>	1 minute	Clean
EN 13697	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i> <i>Escherichia coli</i>	30 seconds	Dirty
EN 16615	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i>	1 minute	Dirty
EN 16615	Bactericidal	<i>Pseudomonas aeruginosa</i> <i>Staphylococcus aureus</i> <i>Enterococcus hirae</i>	1 minute	Clean
EN 14348	Mycobactericidal	<i>Mycobacterium terrae</i> <i>Mycobacterium avium</i>	1 minute	Dirty
EN 14348	Mycobactericidal	<i>Mycobacterium terrae</i> <i>Mycobacterium avium</i>	1 minute	Clean
EN 1650	Fungicidal Yeast (C.a) Fungi (A.b)	<i>Candida albicans</i> <i>Aspergillus brasiliensis</i>	1 minute	Dirty
EN 13624	Yeasticidal	<i>Candida albicans</i>	3 minutes	Clean
EN 13697	Yeasticidal	<i>Candida albicans</i>	1 minute	Clean
EN 16615	Yeasticidal	<i>Candida albicans</i>	3 minutes	Clean

Products within this range



Canister - 200 Wipes

Product Code – W290230

Wipe colour – blue

Packs per case – 10



Bucket - 1000 Wipes

Product Code – W295230

Wipe colour – blue

Packs per case – 2



Bucket - 500 Wipes

Product Code – W84230T

Wipe colour – blue

Packs per case – 2

Manufacturer:
Pal International Limited

Unit 3 Mountpark, Bardon II,
Victoria Road,
Ellistown,
Coalville,
LE67 1FA

T: +44 (0) 1455 555700

E: sales@palinternational.com

W: www.palinternational.com

EU Authorisation Holder:
Pal Hygiene
Products Limited

Unit 5B & Unit 5H,
Fingal Bay Business Park,
Balbriggan, Co. Dublin.
Republic of Ireland

T: +353 15060172

E: sales@palinternational.com

